Membranventil

Metall, DN 4 - 100

Diaphragm Valve

Metal, DN 4 - 100

- Ⓓ ORIGINAL EINBAU- UND MONTAGEANLEITUNG
- ⒼB INSTALLATION, OPERATING AND MAINTENANCE INSTRUCTIONS

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Contents

1	General information	27
2	General safety information	27
2.1	Information for service and operating personnel	28
2.2	Warning notes	28
2.3	Symbols used	29
3	Definition of terms	29
4	Intended area of use	29
5	Technical data	29
6	Order data (2/2-way valves)	33
7	Manufacturer's information	35
7.1	Transport	35
7.2	Delivery and performance	35
7.3	Storage	35
7.4	Tools required	35
8	Functional description	35
9	Construction	35
10	Installation and operation	36
10.1	Installing the diaphragm valve	36
10.2	Control functions	37
10.3	Connecting the control medium	38
10.4	Optical position indicator	38
11	Assembly / disassembly of spare parts	38
11.1	Valve disassembly (removing actuator from body)	38
11.2	Removing the diaphragm	39
11.3	Mounting the diaphragm	39
11.3.1	General information	39
11.3.2	Mounting a concave diaphragm	41
11.3.3	Mounting a convex diaphragm	41
11.4	Actuator mounting on valve body	42
12	Commissioning	42
13	Inspection and servicing	43
14	Disassembly	43
15	Disposal	43
16	Returns	43
17	Information	44
18	Troubleshooting / Fault clearance	45
19	Sectional drawings and spare parts	46
20	Declaration of incorporation	50
21	EU Declaration of conformity	51

1 General information

Prerequisites to ensure that the GEMÜ valve functions correctly:

- x Correct transport and storage
- x Installation and commissioning by trained personnel
- x Operation according to these installation, operating and maintenance instructions
- x Recommended maintenance

Correct installation, operation, servicing and repair work ensure faultless diaphragm valve operation.

The descriptions and instructions apply to the standard versions. For special versions not described in these installation, operating and maintenance instructions the basic information contained herein applies in combination with any additional special documentation.

All rights including copyright and industrial property rights are expressly reserved.

2 General safety information

The safety information does not take into account:

- x Coincidences and events, which may occur during installation, operation and servicing.
- x Local safety regulations which must be adhered to by the operator and by any additional installation personnel.

2.1 Information for service and operating personnel

The installation, operating and maintenance instructions contain fundamental safety information that must be observed during commissioning, operation and servicing. Non-compliance with these instructions may cause:

- x Personal hazard due to electrical, mechanical and chemical effects.
- x Hazard to nearby equipment.
- x Failure of important functions.
- x Hazard to the environment due to the leakage of dangerous materials.

Prior to commissioning:

- Read the installation, operating and maintenance instructions.
- Provide adequate training for the installation and operating personnel.
- Ensure that the contents of the installation, operating and maintenance instructions have been fully understood by the responsible personnel.
- Define the areas of responsibility.

During operation:

- Keep the installation, operating and maintenance instructions available at the place of use.
- Observe the safety information.
- Use only in accordance with the specifications.
- Any servicing work and repairs not described in the installation, operating and maintenance instructions must not be performed without consulting the manufacturer first.

⚠ DANGER

Strictly observe the safety data sheets or the safety regulations that are valid for the media used.

In cases of uncertainty:

- x Consult the nearest GEMÜ sales office.

2.2 Warning notes

Wherever possible, warning notes are organised according to the following scheme:

⚠ SIGNAL WORD

Type and source of the danger

- Possible consequences of non-observance.
- Measures for avoiding danger.

Warning notes are always marked with a signal word and sometimes also with a symbol for the specific danger.

The following signal words and danger levels are used:

⚠ DANGER

Imminent danger!

- Non-observance will lead to death or severe injury.

⚠ WARNING

Potentially dangerous situation!

- Non-observance can cause death or severe injury.

⚠ CAUTION

Potentially dangerous situation!

- Non-observance can cause moderate to light injury.

CAUTION (WITHOUT SYMBOL)

Potentially dangerous situation!

- Non-observance can cause damage to property.

2.3 Symbols used

	Danger - hot surfaces!
	Danger - corrosive materials!
	Hand: indicates general information and recommendations.
	Bullet point: indicates the tasks to be performed.
	Arrow: indicates the response(s) to tasks.
	Enumeration sign

3 Definition of terms

Working medium

The medium that flows through the diaphragm valve.

Control medium

The medium whose increasing or decreasing pressure causes the diaphragm valve to be actuated and operated.

Control function

The possible actuation functions of the diaphragm valve.

5 Technical data

Working medium

Corrosive, inert, gaseous and liquid media which have no negative impact on the physical and chemical properties of the body and diaphragm material.

The valve will seal in both flow directions up to full operating pressure (gauge pressure).

Temperatures

Medium temperature

-10 to 100 °C

Sterilisation temperature ⁽¹⁾

EPDM (code 13/3A)	max. 150 °C ⁽²⁾ , max. 60 min per cycle
EPDM (code 17)	max. 150 °C ⁽²⁾ , max. 180 min per cycle
PTFE/EPDM (code 54)	max. 150 °C ⁽²⁾ , no time limit per cycle
PTFE/EPDM (code 5M)	max. 150 °C ⁽²⁾ , no time limit per cycle

¹ The sterilisation temperature is valid for steam (saturated steam) or superheated water.


² If the sterilisation temperatures listed above are applied to the EPDM diaphragms for longer periods of time, the service life of the diaphragms will be reduced. In these cases, maintenance cycles must be adapted accordingly. This also applies to PTFE diaphragms exposed to high temperature fluctuations.

PTFE diaphragms can also be used as moisture barriers; however, this will reduce their service life.

The maintenance cycles must be adapted accordingly.

GEMÜ 555 and 505 globe valves are particularly suitable for use in the area of steam generation and distribution.

The following valve arrangement for interfaces between steam pipes and process pipes has proven itself over time: A globe valve for shutting off steam pipes and a diaphragm valve as an interface to the process pipes.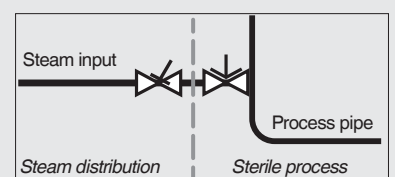

Ambient temperature

0 ... 60 °C

4 Intended area of use

- x The GEMÜ 650 diaphragm valve is designed for installation in piping systems. It controls a flowing medium by being opened or closed by a control medium.
- x **The valve may only be used providing the product technical criteria are complied with (see chapter 5 "Technical data").**
- x Do not paint the bolts and plastic parts of the diaphragm valve!

⚠ WARNING

Use the diaphragm valve only for the intended purpose!

- Otherwise the manufacturer liability and guarantee will be void.
- Use the diaphragm valve only in accordance with the operating conditions specified in the contract documentation and in the installation, operating and maintenance instructions.
- The diaphragm valve may only be used in potentially explosive zones confirmed in the declaration of conformity (ATEX).

Control medium							
Inert gases							
Max. permissible temperature of control medium							60 °C
Filling volume							
Diaphragm size	DN	Actuator size	Actuator version	Spring set	Control function 1	Control function 2	
8	4 to 15	0	T/R	1	0.01 dm ³	0.01 dm ³	
			T/R	A	0.02 dm ³	0.01 dm ³	
10	10 to 20	1	T/R/D/B	1	0.03 dm ³	0.07 dm ³	
25	15 to 25	2	T/R/D/B	1	0.13 dm ³	0.22 dm ³	
40	32 to 40	3	T/R/D/B	1	0.23 dm ³	0.50 dm ³	
			T/R	A	0.50 dm ³	-	
50	50 to 65	4	T/R/D/B	1	0.50 dm ³	1.20 dm ³	
80	65 to 80	5	T/R	1	2.68 dm ³	3.20 dm ³	
			T/R	A	2.13 dm ³	-	
100	100	6	T/R	1	2.78 dm ³	3.40 dm ³	
			T/R	A	2.15 dm ³	-	

C.f. 3 = for filling volume in open position see c.f. 1; for filling volume in closed position see c.f. 2

Operating pressure [bar]								
MG	DN	Control function	Actuator version	EPDM		PTFE		
				Diaphragm material	All valve body materials	Diaphragm material	Forged body	Investment cast body
8	4 to 15	1	0T1, 0R1	3A, 17	0 to 8	5A	0 to 6	0 to 6
			0TA, 0RA		0 to 10		0 to 10	0 to 6
		2 + 3	0T1, 0R1, 0TA, 0RA		0 to 10		0 to 10	0 to 6
10	10 to 20	1	1T1, 1R1	13, 17	0 to 10	54, 5M	0 to 10	0 to 6
			1D1, 1B1		0 to 10		0 to 6	
		2 + 3	1T1, 1R1		0 to 10		0 to 10	0 to 6
			1D1, 1B1		0 to 10		0 to 10	0 to 6
25	15 to 25	1	2T1, 2R1	13, 17	0 to 10	54, 5M	0 to 10	0 to 6
			2D1, 2B1		0 to 10		0 to 6	
		2 + 3	2T1, 2R1		0 to 10		0 to 10	0 to 6
			2D1, 2B1		0 to 10		0 to 10	0 to 6
40	32 to 40	1	3T1, 3R1, 3D1, 3B1	13, 17	0 to 10	54, 5M	0 to 6	0 to 6
			3TA, 3RA		-		0 to 10	0 to 6
		2 + 3	3T1, 3R1		0 to 10		0 to 10	0 to 6
			3D1, 3B1		0 to 10		0 to 10	0 to 6
50	50 to 65	1	4T1, 4R1	13, 17	0 to 10	54, 5M	0 to 10	0 to 6
			4D1, 4B1		0 to 10		0 to 6	
		2 + 3	4T1, 4R1		0 to 10		0 to 10	0 to 6
			4D1, 4B1		0 to 10		0 to 10	0 to 6
80	65 to 80	1	5T1, 5R1	13, 17	0 to 8	54, 5M	0 to 5	-
			5TA, 5RA		-		0 to 10	-
		2 + 3	5T1, 5R1		0 to 8		0 to 5	-
100	100	1	6T1, 6R1	13, 17	0 to 6	54, 5M	0 to 4	-
			6TA, 6RA		-		0 to 10	-
		2 + 3	6T1, 6R1		0 to 6		0 to 4	-

All pressures are gauge pressures. Operating pressure values were determined with static operating pressure applied on one side of a closed valve. Sealing at the valve seat and atmospheric sealing is ensured for the given values.
Information on operating pressures applied on both sides and for high purity media on request
MG = diaphragm size

Control pressure [bar]				
MG	DN	Control function	Actuator version	Control pressure
8	4 to 15	1	0T1, 0R1	5,0 to 7,0
			0TA, 0RA	3,5 to 7,0
		2 + 3	0T1, 0R1	max. 5,5
			0TA, 0RA	max. 4,5
10	10 to 20	1	1T1, 1R1, 1D1, 1B1	4,5 to 7,0
		2 + 3	1T1, 1R1, 1D1, 1B1	max. 4,5
25	15 to 25	1	2T1, 2R1, 2D1, 2B1	5,0 to 7,0
		2 + 3	2T1, 2R1, 2D1, 2B1	max. 4,5
40	32 to 40	1	3T1, 3R1, 3D1, 3B1	4,5 to 7,0
			3TA, 3RA	3,5 to 7,0
		2 + 3	3T1, 3R1, 3D1, 3B1	max. 5,5
50	50 to 65	1	4T1, 4R1, 4D1, 4B1	4,5 to 7,0
		2 + 3	4T1, 4R1, 4D1, 4B1	max. 4,5
80	65 to 80	1	5T1, 5R1	3,5 to 7,0
			5TA, 5RA	4,5 to 7,0
		2 + 3	5T1, 5R1	max. 3,5
100	100	1	6T1, 6R1	3,5 to 7,0
			6TA, 6RA	5,0 to 7,0
		2 + 3	6T1, 6R1	max. 3,5

MG = diaphragm size

Kv values [m ³ /h]									
Pipe standard		DIN	EN 10357 series B (formerly DIN 11850 series 1)	EN 10357 series A (formerly DIN 11850 series 2) / DIN 11866 series A	DIN 11850 Series 3	SMS 3008	ASME BPE / DIN 11866 series C	ISO 1127 / EN 10357 series C / DIN 11866 series B	DIN ISO 228
Connection code		0	16	17	18	37	59	60	1
MG	DN								
8	4	0.5	-	-	-	-	-	-	-
	6	-	-	1.1	-	-	-	1.2	-
	8	-	-	1.3	-	-	0.6	2.2	-
	10	-	2.1	2.1	2.1	-	1.3	-	-
	15	-	-	-	-	-	2.0	-	-
10	10	-	2.4	2.4	2.4	-	2.2	3.3	-
	12	-	-	-	-	-	-	-	3.2
	15	3.3	3.8	3.8	3.8	-	2.2	4.0	3.4
	20	-	-	-	-	-	3.8	-	-
25	15	4.1	4.7	4.7	4.7	-	-	7.4	6.5
	20	6.3	7.0	7.0	7.0	-	4.4	13.2	10.0
	25	13.9	15.0	15.0	15.0	12.6	12.2	16.2	14.0
40	32	25.3	27.0	27.0	27.0	26.2	-	30.0	26.0
	40	29.3	30.9	30.9	30.9	30.2	29.5	32.8	33.0
50	50	46.5	48.4	48.4	48.4	51.7	50.6	55.2	60.0
	65	-	-	-	-	62.2	61.8	-	-
80	65	-	-	77.0	-	68.5	68.5	96.0	-
	80	-	-	111.0	-	80.0	87.0	111.0	-
100	100	-	-	194.0	-	173.0	188.0	214.0	-

MG = diaphragm size


Kv values determined acc. to DIN EN 60534, inlet pressure 5 bar, Δp 1 bar, stainless steel valve body (forged body) and soft elastomer diaphragm. The Kv values for other product configurations (e.g. other diaphragm or body materials) may differ. In general, all diaphragms are subject to the influences of pressure, temperature, the process and their tightening torques. Therefore the Kv values may exceed the tolerance limits of the standard.

The Kv value curve (Kv value dependent on valve stroke) can vary depending on the diaphragm material and duration of use.


Autoclavability

Actuator size 0	Standard version with autoclave capability
Actuator size 1	Standard version with autoclave capability
Actuator size 2	Standard version with autoclave capability
Actuator size 3	with special version
Actuator size 4	with special version
Actuator size 5	not possible
Actuator size 6	not possible


**Control function 2 + 3
with elastomer diaphragm
Diaphragm size 8 - 25**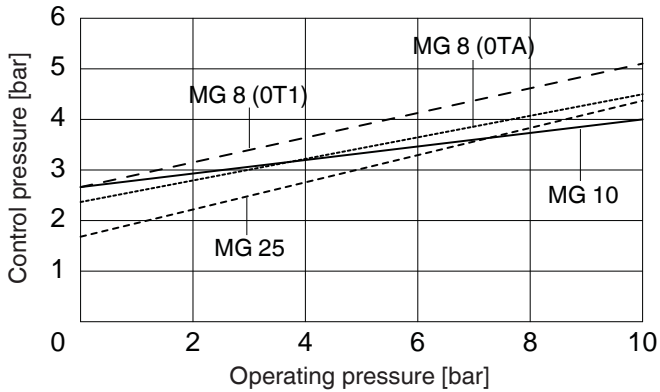


**Control function 2 + 3
with elastomer diaphragm
Diaphragm size 40 - 100**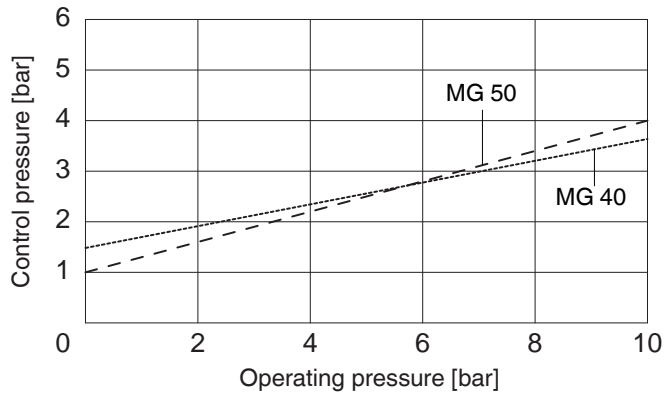

**Control function 2 + 3
with PTFE diaphragm
Diaphragm size 8 - 25**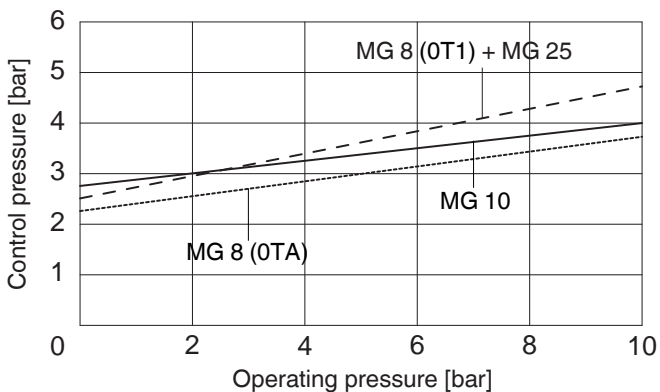

**Control function 2 + 3
with PTFE diaphragm
Diaphragm size 40 - 100**

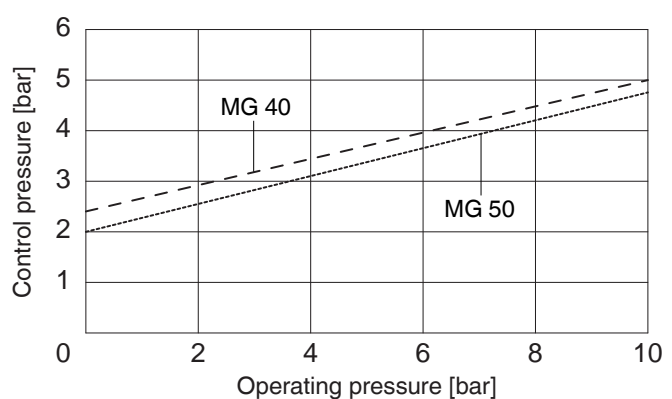The control pressure depending on the prevailing operating pressure, as shown in the diagram, is intended as a guide for operating the system with low wear on the diaphragm.

6 Order data (2/2-way valves)

Body configuration	Code
Tank bottom valve body (actuator version T)	B**
2/2-way body (actuator version D and T)	D
T body (actuator version T)	T*
* For dimensions see T Valves brochure	
** Dimensions and versions on request	

Connection	Code
Butt weld spigots	
Spigots DIN	0
Spigots EN 10357 series B (formerly DIN 11850 series 1)	16
Spigot EN 10357 series A (formerly DIN 11850 series 2) / DIN 11866 series A	17
Spigots DIN 11850 series 3	18
Spigots JIS-G 3447	35
Spigots JIS-G 3459	36
Spigots SMS 3008	37
Spigots BS 4825 Part 1	55
Spigot ASME BPE / DIN 11866 series C	59
Spigot ISO 1127 / EN 10357 series C / DIN 11866 series B	60
Spigots ANSI/ASME B36.19M Schedule 10s	63
Spigots ANSI/ASME B36.19M Schedule 40s	65
Threaded connections	
Threaded sockets DIN ISO 228	1
Threaded spigots DIN 11851	6
One side threaded spigot, other side cone spigot and union nut, DIN 11851	62
Aseptic unions on request	
Flanges	
Flanges EN 1092 / PN16 / form B, length EN 558, series 1, ISO 5752, basic series 1	8*
Flanges ANSI Class 150 RF, length MSS SP-88	38*
Flanges ANSI Class 125/150 RF, length EN 558, series 1, ISO 5752, basic series 1	39*
Clamp connections	
Clamps ASME BPE for pipe ASME BPE, length ASME BPE	80
Clamps DIN 32676 series B for pipe EN ISO 1127, length EN 558, series 7	82
Clamps ASME BPE for pipe ASME BPE, length EN 558, series 7	88
Clamps DIN 32676 series A for pipe DIN 11850, length EN 558, series 7	8A
Clamps SMS 3017 for pipe SMS 3008, length EN 558, series 7	8E
Aseptic clamps on request	
* Connection code 8, 38, 39 only possible in conjunction with actuator version code B / R	
For overview of available valve bodies see page 15/16	

Valve body material	Code
1.4435, investment casting	C3
1.4408, investment casting	37
1.4408, PFA lined	39
1.4435 (316L), forged body	40
1.4435 (BN2), forged body Δ Fe<0.5%	42
1.4539, forged body	F4

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Control function	Code
Normally closed (NC)	1
Normally open (NO)	2
Double acting (DA) (with opening spring)	3

Actuator size	Code
Actuator size 0 (diaphragm size 8)	0
Actuator size 1 (diaphragm size 10)	1
Actuator size 2 (diaphragm size 25)	2
Actuator size 3 (diaphragm size 40)	3
Actuator size 4 (diaphragm size 50)	4
Actuator size 5 (diaphragm size 80)	5
Actuator size 6 (diaphragm size 100)	6

Actuator version	Code
Only for body configuration D (diaphragm size 10 - 50)	D
Only for body configuration D (diaphragm size 10 - 50) Control medium connector 90° offset to flow direction	B
For body config. B, D, M and T (diaphragm size 8 - 100)	T
For body config. B, D, M and T (diaphragm size 8 - 100) Control medium connector 90° offset to flow direction	R

Spring set	Code
Standard	1
For higher operating pressure (diaphragm size 8)	A

Internal surface finishes for forged and block material bodies ¹

Readings for Process Contact Surfaces	Mechanically polished ²		Electropolished	
	Hygienic class DIN 11866	Code	Hygienic class DIN 11866	Code
Ra ≤ 0.80 µm	H3	1502	HE3	1503
Ra ≤ 0.60 µm	-	1507	-	1508
Ra ≤ 0.40 µm	H4	1536	HE4	1537
Ra ≤ 0.25 µm ³	H5	1527	HE5	1516

Readings for Process Contact Surfaces acc. to ASME BPE 2016 ⁴	Mechanically polished ²		Electropolished	
	ASME BPE Surface Designation	Code	ASME BPE Surface Designation	Code
Ra Max. = 0.76 µm (30 µinch)	SF3	SF3	-	-
Ra Max. = 0.64 µm (25 µinch)	SF2	SF2	SF6	SF6
Ra Max. = 0.51 µm (20 µinch)	SF1	SF1	SF5	SF5
Ra Max. = 0.38 µm (15 µinch)	-	-	SF4	SF4

Internal surface finishes for investment cast bodies

Readings for Process Contact Surfaces	Mechanically polished ²	
	Hygienic class DIN 11866	Code
Ra ≤ 6.30 µm	-	1500
Ra ≤ 0.80 µm	H3	1502
Ra ≤ 0.60 µm ⁵	-	1507

¹ Surface finishes of customized valve bodies may be limited in special cases.

² Or any other finishing method that meets the Ra value (acc. to ASME BPE).

³ The smallest possible Ra finish for pipe connections with an internal pipe diameter < 6 mm is 0.38 µm.

⁴ When using these surfaces, the bodies are marked according to the specifications of ASME BPE.

The surfaces are only available for valve bodies which are made of materials (e.g. GEMÜ material codes 40, 41, F4, 44) and use connections (e.g. GEMÜ connection codes 59, 80, 88) according to ASME BPE.

⁵ Not possible for GEMÜ connection code 59, DN 8 and GEMÜ connection code 0, DN 4.

Ra acc. to DIN EN ISO 4288 and ASME B46.1

Special function	Code
3-A compliant design	M

Order example	650	50	D	60	40	54	1	4	T	1	1503	M
Type	650											
Nominal size		50										
Body configuration (code)			D									
Connection (code)				60								
Valve body material (code)					40							
Diaphragm material (code)						54						
Control function (code)							1					
Actuator size (code)								4				
Actuator version (code)									T			
Spring set (code)										1		
Surface finish (code)											1503	
Special function (Code)												M

7 Manufacturer's information

7.1 Transport

- Only transport the diaphragm valve by suitable means. Do not drop. Handle carefully.
- Dispose of packing material according to relevant local or national disposal regulations / environmental protection laws.

7.2 Delivery and performance

- Check that all parts are present and check for any damage immediately upon receipt.
- The scope of delivery is apparent from the dispatch documents and the design from the order number.
- The valve's delivery condition:

Control function:	Condition:
1 Normally closed (NC)	closed
2 Normally open (NO)	open
3 Double acting (DA)	open

- The performance of the diaphragm valve is checked at the factory.

7.3 Storage

- Store the diaphragm valve free from dust and moisture in its original packaging.
- Avoid UV rays and direct sunlight.
- Maximum storage temperature: 40 °C.
- Solvents, chemicals, acids, fuels or similar fluids must not be stored in the same room as valves and their spare parts.


7.4 Tools required

- The tools required for installation and assembly are **not** included in the scope of delivery.
- Use appropriate, functional and safe tools.

8 Functional description

GEMÜ 650 is a metal diaphragm valve with a 2/2-way straight through body, T or tank bottom valve body or in multi-port design. The valve has a piston actuator and an optical position indicator as standard. All metallic actuator components are made of stainless steel. The compression springs of diaphragm sizes 80 and 100 are made of epoxy coated spring steel. "Normally closed (NC)", "Normally open (NO)" and "Double acting (DA)" control functions are available. The valve body and the diaphragm are available in various designs as shown in the data sheet. The valve has CIP / SIP cleaning and sterilising capabilities (autoclave capability dependent on version). Diverse accessories are available, such as stroke limiters, electrical position indicators, positioners.

9 Construction

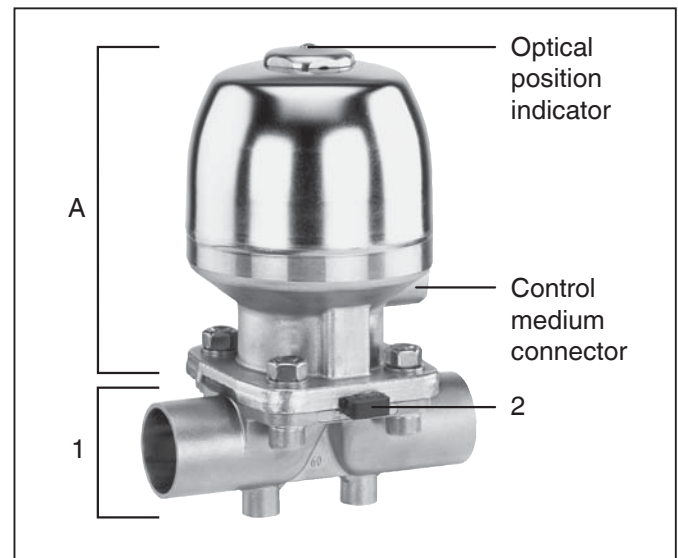Construction

1	Valve body
2	Diaphragm
A	Actuator

10 Installation and operation

Prior to installation:

- Ensure that valve body and diaphragm material are appropriate and compatible to handle the working medium.
- **Check the suitability prior to the installation.**
See chapter 5 "Technical data".

10.1 Installing the diaphragm valve

⚠ WARNING

The equipment is subject to pressure!

- Risk of severe injury or death!
- Only work on depressurized plant.

⚠ WARNING

Corrosive chemicals!

- Risk of caustic burns!
- Wear appropriate protective gear when installing.

⚠ CAUTION

Hot plant components!

- Risk of burns!
- Only work on plant that has cooled down.

⚠ CAUTION

Never use the valve as a step or an aid for climbing!

- This entails the risk of slipping-off or damaging the valve.

CAUTION

Do not exceed the maximum permissible pressure!

- Take precautionary measures to avoid possible pressure surges (water hammer).

- Installation work must only be performed by trained personnel.
- Use appropriate protective gear as specified in plant operator's guidelines.

Installation location:

⚠ CAUTION

- Do not apply external force to the valve.
- Choose the installation location so that the valve cannot be used as a foothold (climbing aid).
- Lay the pipeline so that the valve body is protected against transverse and bending forces, and also vibrations and tension.
- Only mount the valve between matching aligned pipes.

x Direction of the working medium: optional.

x Mounting position of the diaphragm valve: optional.

Installation:

1. Ensure the suitability of the valve for each respective use. The valve must be appropriate for the piping system operating conditions (medium, medium concentration, temperature and pressure) and the prevailing ambient conditions. Check the technical data of the valve and the materials.
2. Shut off plant or plant component.
3. Secure against recommissioning.
4. Depressurize the plant or plant component.
5. Completely drain the plant (or plant component) and let it cool down until the temperature is below the media vaporization temperature and scalding can be ruled out.
6. Correctly decontaminate, rinse and ventilate the plant or plant component.

Installation - Butt weld spigots:

1. Adhere to good welding practices!
2. Disassemble the actuator with the diaphragm before welding the valve body into the pipeline (see chapter 11.1).
3. Allow butt weld spigots to cool down.
4. Reassemble the valve body and the actuator with diaphragm (see chapter 11.4).

Installation - Clamp connections:

- When installing clamp connections,

insert a gasket between the body clamp and the adjacent piping clamp and join them using the appropriate clamp fitting. The gasket and the clamp for clamp connections are not included in the scope of delivery.


	<p>Important: Butt weld spigots / Clamp connections: Angle of rotation for welding into pipeline to enable optimised draining see brochure "Angle of rotation for 2/2-way valve bodies" (on request or www.gemu-group.com).</p>
---	--

Installation - Threaded connections:

- Screw the threaded connections into the piping in accordance with valid standards.
- Screw the diaphragm valve body into the piping, use appropriate thread sealant. The thread sealant is not included in the scope of delivery.

Installation - Flange connection:

1. Pay attention to clean, undamaged sealing surfaces on the mating flanges.
2. Align flanges carefully before installing them.
3. Centre the seals accurately.
4. Connect the valve flange and the piping flange using appropriate sealing material and matching bolting. Sealing material and bolts are not included in the scope of delivery.
5. Use all flange holes.
6. Only use connector elements made of approved materials!
7. Tighten the bolts diagonally!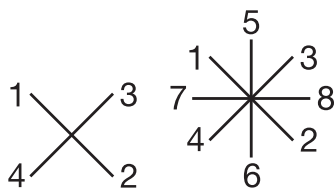

Observe appropriate regulations for connections!

After the installation:

- Reactivate all safety and protective devices.

10.2 Control functions

The following control functions are available:

Control function 1

Normally closed (NC):

Valve resting position: closed by spring force. Activation of the actuator (connector 2) opens the valve. When the actuator is vented, the valve is closed by spring force.

Control function 2



Normally open (NO):

Valve resting position: opened by spring force. Activation of the actuator (connector 4) closes the valve. When the actuator is vented, the valve is opened by spring force.

Control function 3

Double acting (DA):

Valve resting position: opened by spring force. The valve is opened and closed by activating the respective control medium connectors (connector 2: open / connector 4: close).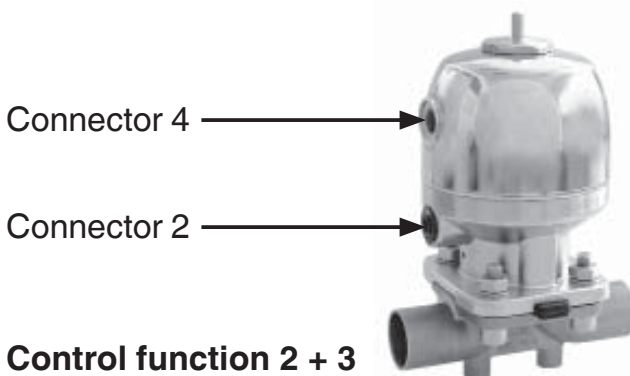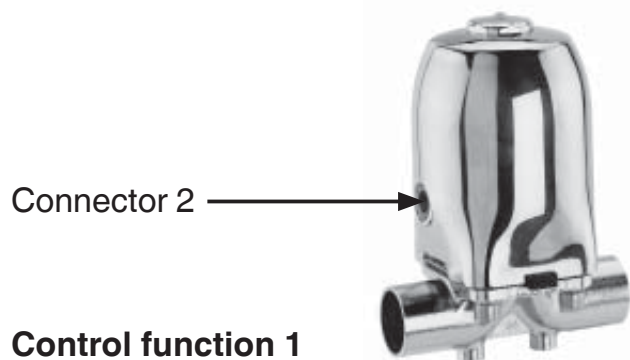

Control function	Connectors	
	2	4
1 (NC)	+	-
2 (NO)	-	+
3 (DA)	+	+
+ = available / - = not available (for connectors 2 / 4 see pictures above)		

10.3 Connecting the control medium

 **Important:**
Connect the control medium lines tension-free and without any bends or knots!
Use appropriate connectors according to the application.


Thread size of the control medium connectors:

Diaphragm size 8: G1/8


Diaphragm size 10 - 100: G1/4

Control function		Connectors
1	Normally closed (NC)	2: Control medium (open)
2	Normally open (NO)	4: Control medium (close)
3	Double acting (DA)	2: Control medium (open) 4: Control medium (close)
For connectors 2 / 4 see pictures on page 34		

10.4 Optical position indicator

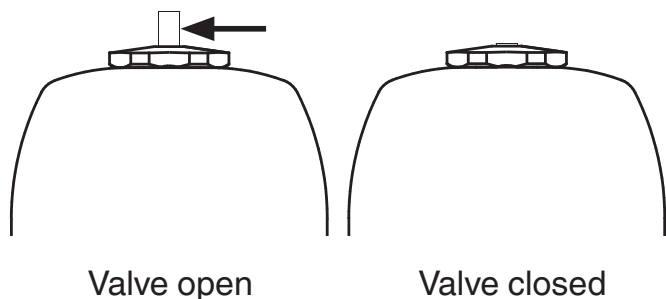
11 Assembly / disassembly of spare parts

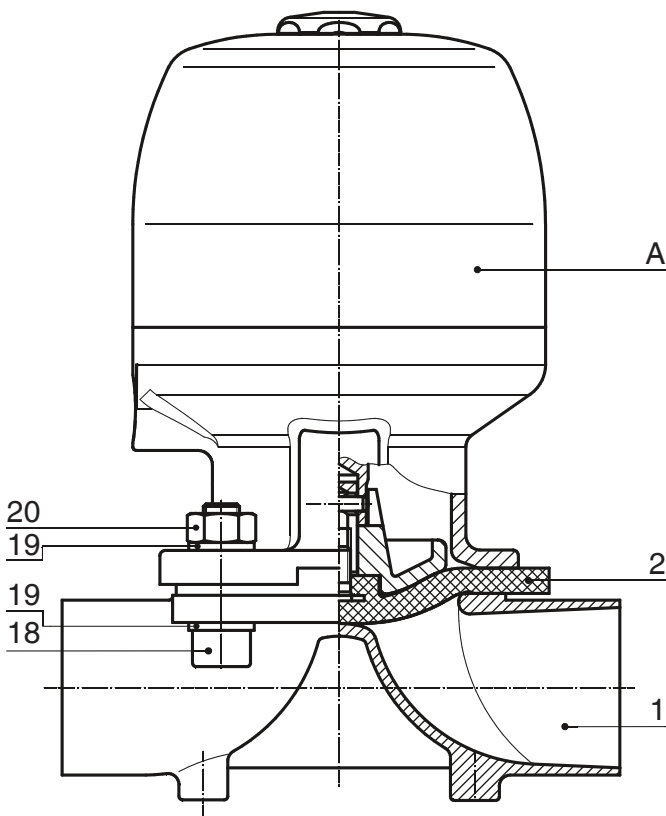11.1 Valve disassembly (removing actuator from body)

1. Move actuator **A** to the open position.
2. Diagonally loosen and remove the fastening elements between valve body **1** and actuator **A**.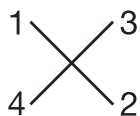

3. Lift off actuator **A** from valve body **1**.
4. Move actuator **A** to the closed position.

 **Important:**
After disassembly, clean all parts of contamination (do not damage parts). Check parts for potential damage, replace if necessary (only use genuine parts from GEMÜ).

11.2 Removing the diaphragm

	Important: Before removing the diaphragm, please remove the actuator, see "Valve disassembly (removing actuator from body)".
---	--

1. Unscrew the diaphragm or pull it out (diaphragm size 8).
2. Clean all parts of the remains of product and contamination. Do not scratch or damage parts during cleaning!
3. Check all parts for potential damage.
4. Replace damaged parts (only use genuine parts from GEMÜ).

11.3 Mounting the diaphragm

11.3.1 General information

	Important: Mount the correct diaphragm that suits the valve (suitable for medium, medium concentration, temperature and pressure). The diaphragm is a wearing part. Check the technical condition and function of the diaphragm valve before commissioning and during the whole term of use. Carry out checks regularly and determine the check intervals in accordance with the conditions of use and / or the regulatory codes and provisions applicable for this application.
--	--

	Important: If the diaphragm is not screwed into the adapter far enough, the closing force is transmitted directly onto the diaphragm pin and not via the compressor. This will cause damage and early failure of the diaphragm and thus leakage of the valve. If the diaphragm is screwed in too far no perfect sealing at the valve seat will be achieved and the function of the valve is no longer ensured.
--	--

	Important: Incorrectly mounted diaphragm may cause valve leakage / emission of medium. In this case remove the diaphragm, check the complete valve and diaphragm and reassemble again proceeding as described above.
---	--

Diaphragm sizes 10 - 80:


The compressor is loose.

Diaphragm sizes 8 and 100:

The compressor is fixed to the spindle.


Diaphragm size 8:

Compressor and actuator flange seen from below: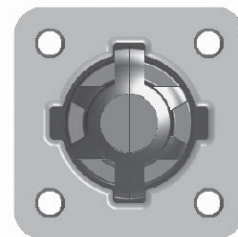

Diaphragm size 10:

Compressor and actuator flange seen from below: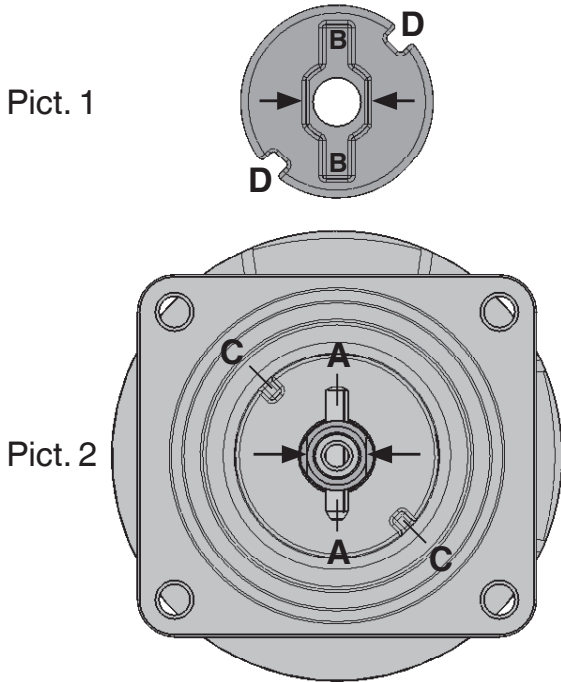

Anti-twist system of the spindle at the compressor


A double flat (arrows picture 2) is fitted at the end of the actuator spindle to protect the spindle against twisting. When mounting the compressor, the double flat must be in correct alignment with the recess of the compressor back (arrows picture 1).

If the actuator spindle is not in the correct position, it must be turned to the correct position. The position of **A** is offset by 45° to the position of **C**.

Place the compressor loosely on the actuator spindle, fit the recesses **D** into the guides **C** and **A** into **B**. The compressor must be able to be moved freely between the guides!

Diaphragm size 25 - 80:


Compressor and actuator flange seen from below: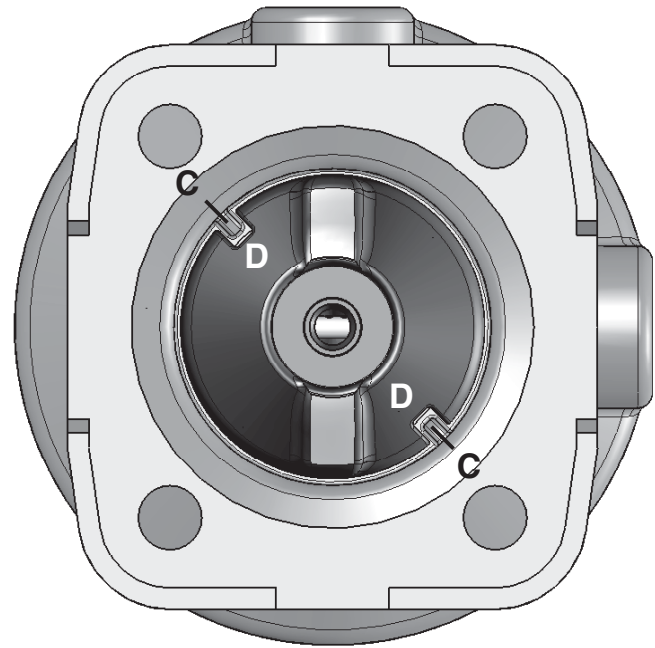

Place the compressor loosely on the actuator spindle, fit the grooves **D** into the guides **C**. The compressor must be able to be moved freely between the guides!

Diaphragm size 100:


Compressor and actuator flange seen from below: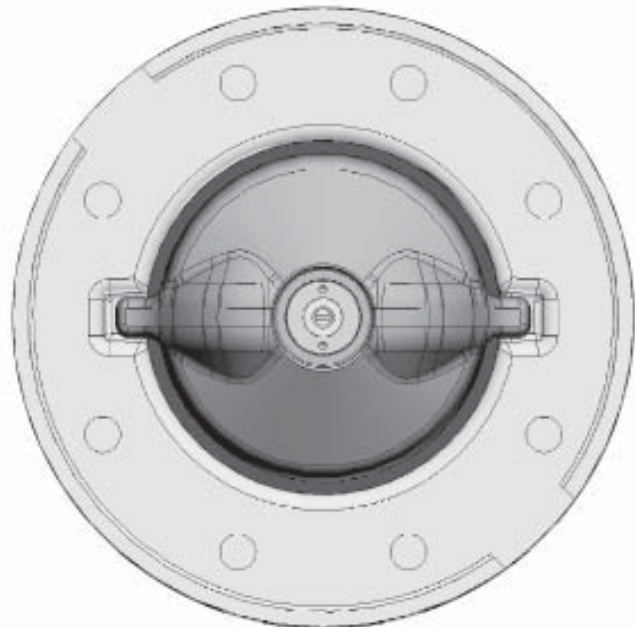

11.3.2 Mounting a concave diaphragm

Diaphragm size 8

Push-fit diaphragm:

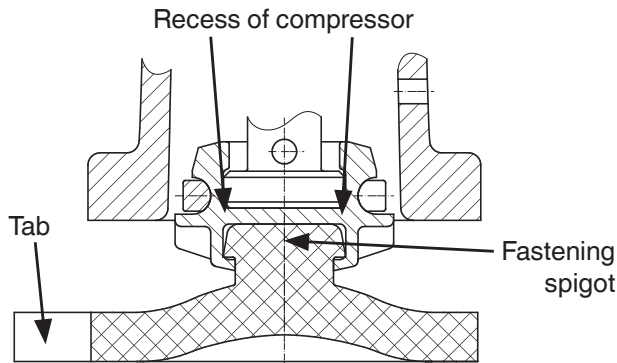1. Move actuator **A** to the closed position.
2. Place the diaphragm **2** with the fastening spigot in an inclined position at the recess of the compressor.


Important:

Do not use any greases or lubricants!

3. Turn the diaphragm as manual force is applied to push the spigot into the compressor.
4. Align diaphragm tab (identifying manufacturer and material) in parallel to compressor weir.

Diaphragm sizes 10 - 100

Threaded pin type diaphragm: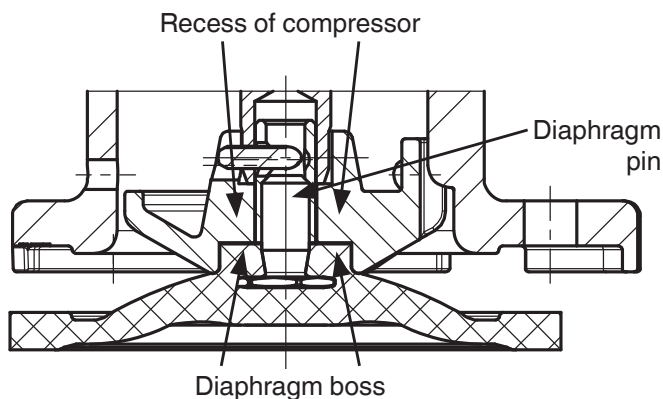


1. Move actuator **A** to the closed position.
2. Diaphragm size 10: ensure that the anti-twist system is engaged. Diaphragm sizes 25 - 80: place the compressor loosely on the actuator spindle, fit the grooves into the guides (see chapter 11.3.1 "General information").


3. Check if the compressor fits closely in the guides.
4. Screw new diaphragm tightly into the compressor manually.
5. Check if the diaphragm boss fits closely in the recess of the compressor.
6. If it is difficult to screw it in, check the thread, replace damaged parts (only use genuine parts from GEMÜ).
7. When clear resistance is felt turn back the diaphragm anticlockwise until its bolt holes are in correct alignment with the bolt holes of the actuator.

11.3.3 Mounting a convex diaphragm

1. Move actuator **A** to the closed position.
2. Diaphragm sizes 25 - 80: place the compressor loosely on the actuator spindle, fit the grooves into the guides (see chapter 11.3.1 "General information").
3. Check if the compressor fits closely in the guides.
4. Invert the new diaphragm face manually; use a clean, padded mat with bigger nominal sizes.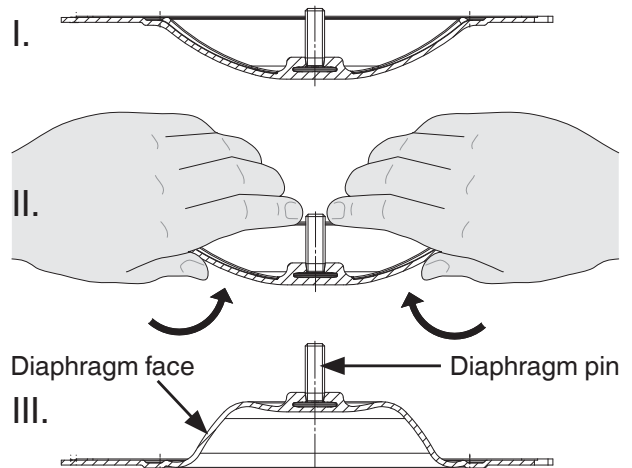

5. Position the new backing diaphragm onto the compressor.
6. Position the diaphragm face onto the backing diaphragm.
7. Screw diaphragm face tightly into the compressor manually. The diaphragm boss must fit closely in the recess of the compressor.

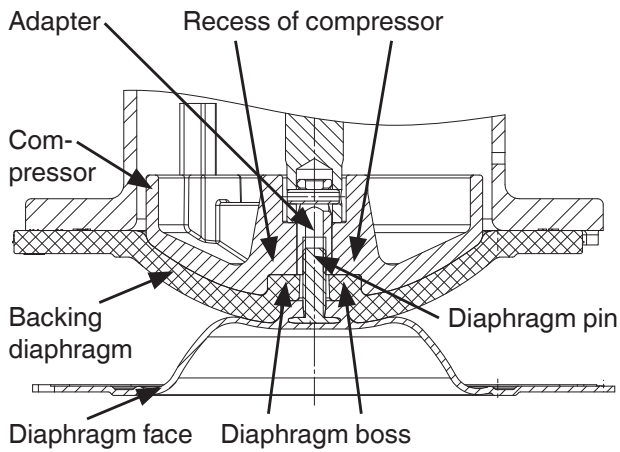8. If it is difficult to screw it in, check the thread, replace damaged parts.
9. When clear resistance is felt turn back the diaphragm anticlockwise until its bolt holes are in correct alignment with the bolt holes of the actuator.
10. Press the diaphragm face tightly onto the backing diaphragm manually so that it returns to its original shape and fits closely on the backing diaphragm.

11.4 Actuator mounting on valve body

1. Move actuator **A** to the open position.
2. Position actuator **A** with the mounted diaphragm **2** on the valve body **1**, take care to align the compressor weir and valve body weir (only for diaphragm size 8).
3. Tighten bolts **18**, washers **19** and nuts **20** by hand (hand tight only) (fastening elements may vary dependent on diaphragm size and / or valve body version).
4. Move actuator **A** to the closed position.
5. Fully tighten the bolts **18** with nuts **20** diagonally.
6. Ensure that the diaphragm **2** is compressed evenly (approx. 10-15 %, visible by an even bulge to the outside).
7. Check tightness of completely assembled valve.

Important:

Service and maintenance: Diaphragms set in the course of time. After valve disassembly / assembly check that the bolts **18** and nuts **20** on the body are tight and retighten as necessary (at the very latest after the first sterilisation process).

12 Commissioning

⚠ WARNING

Corrosive chemicals!

- Risk of caustic burns!
- Check the tightness of the media connections prior to commissioning!
- Use only the appropriate protective gear when performing the tightness check.

⚠ CAUTION

Protect against leakage!

- Provide precautionary measures against exceeding the maximum permitted pressures caused by pressure surges (water hammer).

Prior to cleaning or commissioning the plant:

- Check the tightness and the function of the diaphragm valve (close and reopen the diaphragm valve).
- If the plant is new and after repairs rinse the piping system with a fully opened diaphragm valve (to remove any harmful foreign matter).

Cleaning:

- x The plant operator is responsible for selecting the cleaning material and performing the procedure.

**Important:**

Service and maintenance:
Diaphragms set in the course of time. After valve disassembly / assembly check that the bolts **18** and nuts **20** on the body are tight and retighten as necessary (at the very latest after the first sterilisation process).

13 Inspection and servicing

⚠ WARNING**The equipment is subject to pressure!**

- Risk of severe injury or death!
- Only work on depressurized plant.

⚠ CAUTION**Hot plant components!**

- Risk of burns!
- Only work on plant that has cooled down.

⚠ CAUTION

- Servicing and maintenance work may only be performed by trained personnel.
- GEMÜ shall assume no liability whatsoever for damages caused by improper handling or third-party actions.
- In case of doubt, contact GEMÜ before commissioning.

1. Use appropriate protective gear as specified in plant operator's guidelines.
2. Shut off plant or plant component.
3. Secure against recommissioning.
4. Depressurize the plant or plant component.

The operator must carry out regular visual examination of the valves dependent on the operating conditions and the potential danger in order to prevent leakage and damage. The valve also has to be disassembled in the corresponding intervals and checked for wear (see chapter 11 "Assembly / Disassembly of spare parts").

14 Disassembly

Disassembly is performed observing the same precautionary measures as for installation.

- Disassemble the diaphragm valve (see chapter 11.1 "Valve disassembly (removing actuator from body)").

15 Disposal

- All valve parts must be disposed of according to relevant local or national disposal regulations / environmental protection laws.
- Pay attention to adhered residual material and gas diffusion from penetrated media.

16 Returns

- Clean the diaphragm valve.
- Request a goods return declaration form from GEMÜ.
- Returns must be made with a completed declaration of return.

If not completed, GEMÜ cannot process
x credits or
x repair work
but will dispose of the goods at the operator's expense.

**Note for returns:**

Legal regulations for the protection of the environment and personnel require that you include the completed and signed goods return declaration with the dispatch documents. Your returned goods can be processed only when this declaration is completed.

17 Information

	<p>Note on Directive 2014/34/EU (ATEX Directive): A supplement to Directive 2014/34/EU is included with the product if it was ordered according to ATEX.</p>
	<p>Note on staff training: Please contact us at the address on the last page for staff training information.</p>


Should there be any doubts or misunderstandings in the preceding text, the German version of this document is the authoritative document!

18 Troubleshooting / Fault clearance


Fault	Possible cause	Fault clearance
Control medium escapes from vent hole* / vent* in the actuator cover (control function NC) or from connector 2 (control function NO (see chapter 10.2 "Control functions"))	Actuator piston faulty	Replace actuator
Control medium escapes from leak detection hole*	Spindle seal leaking	Replace actuator and check control medium for impurities
Working medium escapes from leak detection hole*	Valve diaphragm faulty	Check valve diaphragm for damage, replace diaphragm if necessary
Valve doesn't open or doesn't open fully	Control pressure too low (for control function NC)	Operate valve with control pressure specified in data sheet
	Pilot valve faulty	Check and replace pilot valve
	Actuator faulty	Replace actuator
	Control medium not connected	Connect control medium
	Valve diaphragm incorrectly mounted	Remove actuator, check diaphragm mounting, replace if necessary
	Actuator spring faulty (for control function NO)	Replace actuator
Valve leaks downstream (doesn't close or doesn't close fully)	Operating pressure too high	Operate valve with operating pressure specified in data sheet
	Control pressure too low (for control function NO and control function DA)	Operate valve with control pressure specified in data sheet
	Foreign matter between valve diaphragm and valve body weir	Remove actuator, remove foreign matter, check valve diaphragm and valve body weir for damage and replace if necessary
	Valve body weir leaking or damaged	Check valve body weir for damage, if necessary replace valve body
	Valve diaphragm faulty	Check valve diaphragm for damage, replace diaphragm if necessary
	Actuator spring faulty (for control function NC)	Replace actuator
Valve leaks between actuator and valve body	Valve diaphragm incorrectly mounted	Remove actuator, check diaphragm mounting, replace if necessary
	Bolting between valve body and actuator loose	Retighten bolting between valve body and actuator
	Valve diaphragm faulty	Check valve diaphragm for damage, replace diaphragm if necessary
	Actuator / valve body damaged	Replace actuator / valve body
Valve body connection to piping leaks	Incorrect installation	Check installation of valve body in piping
	Bolting / threaded connections loose	Tighten bolting / threaded connections
	Sealing material faulty	Replace sealing material
Valve body leaks	Valve body faulty or corroded	Check valve body for damage, replace valve body if necessary

* see chapter 19 "Sectional drawings and spare parts"


19 Sectional drawings and spare parts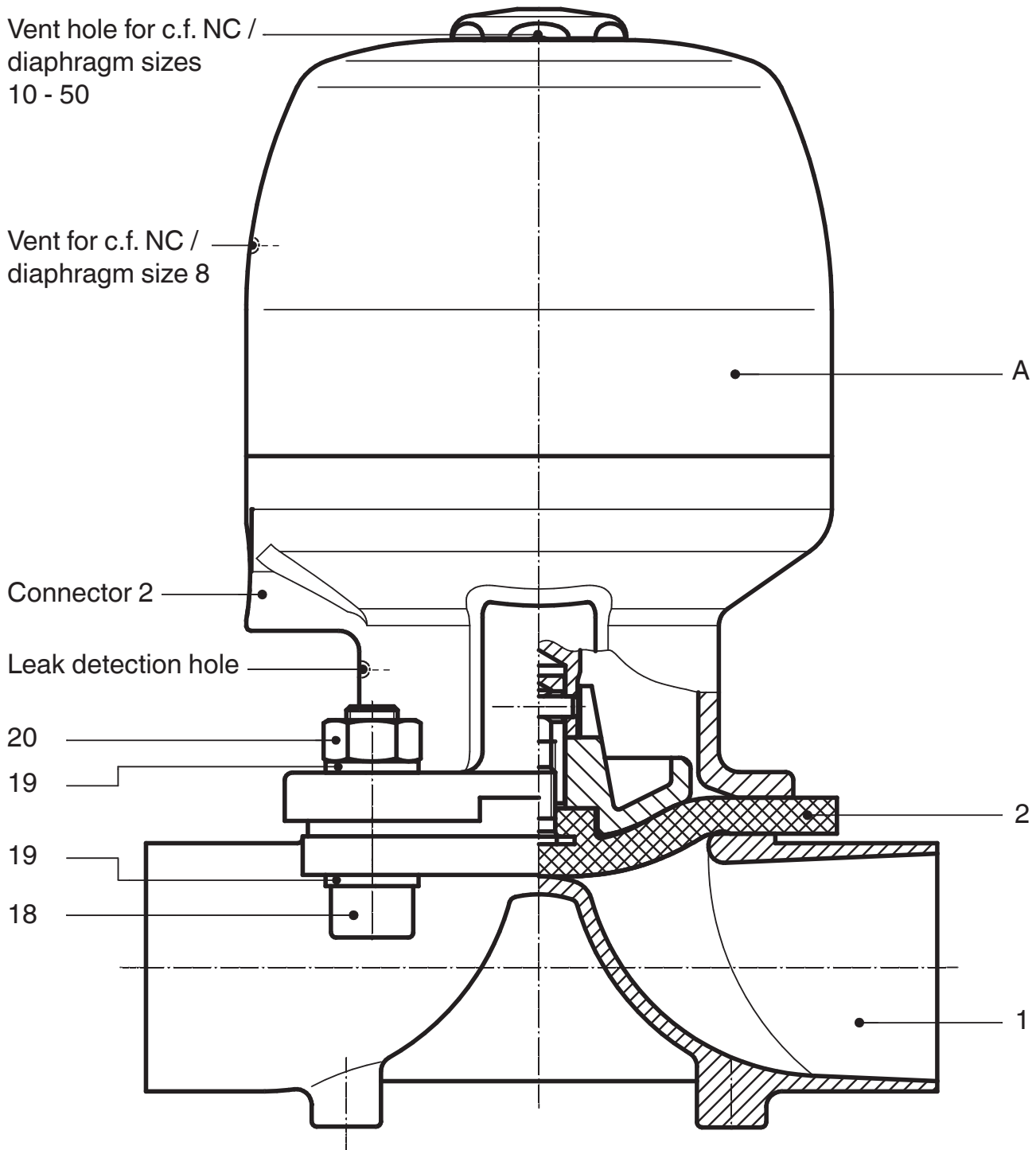


Item	Name	Order description
Diaphragm sizes 8 - 50 / Actuator version: T		
1	Valve body	K600...
2	Diaphragm	600...M
18	Bolt	} 650...S30...
19	Washer	
20	Nut	
A	Actuator	9650...

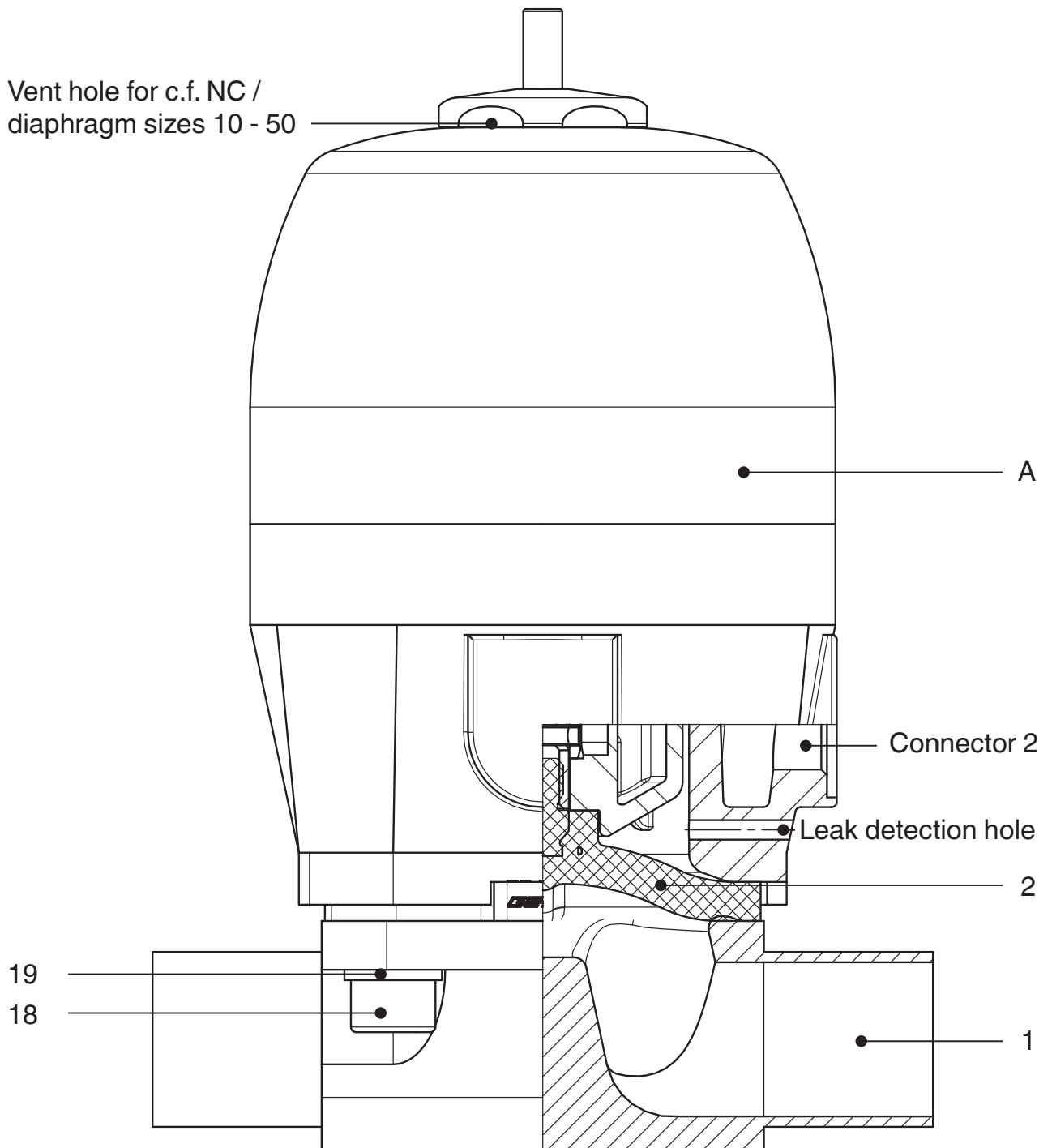Item	Name	Order description
Diaphragm sizes 10 - 50 / Actuator version: D		
1	Valve body	K600...
2	Diaphragm	600...M
18	Bolt	} 650...S30...
19	Washer	
A	Actuator	9650...

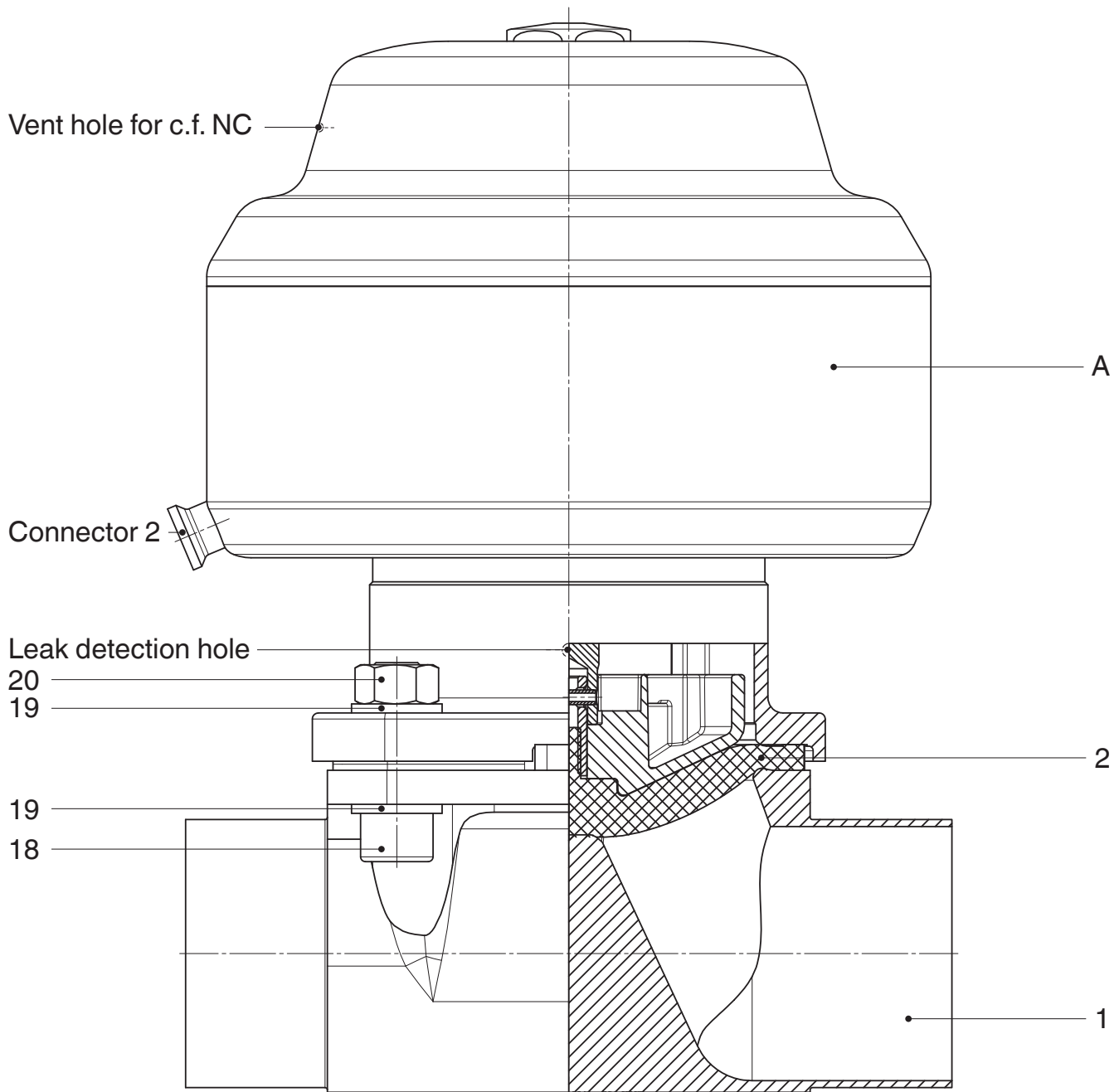Item	Name	Order description
Diaphragm size 80		
1	Valve body	K600...
2	Diaphragm	600...M
18	Bolt	} 650...S30...
19	Washer	
20	Nut	
A	Actuator	9650...

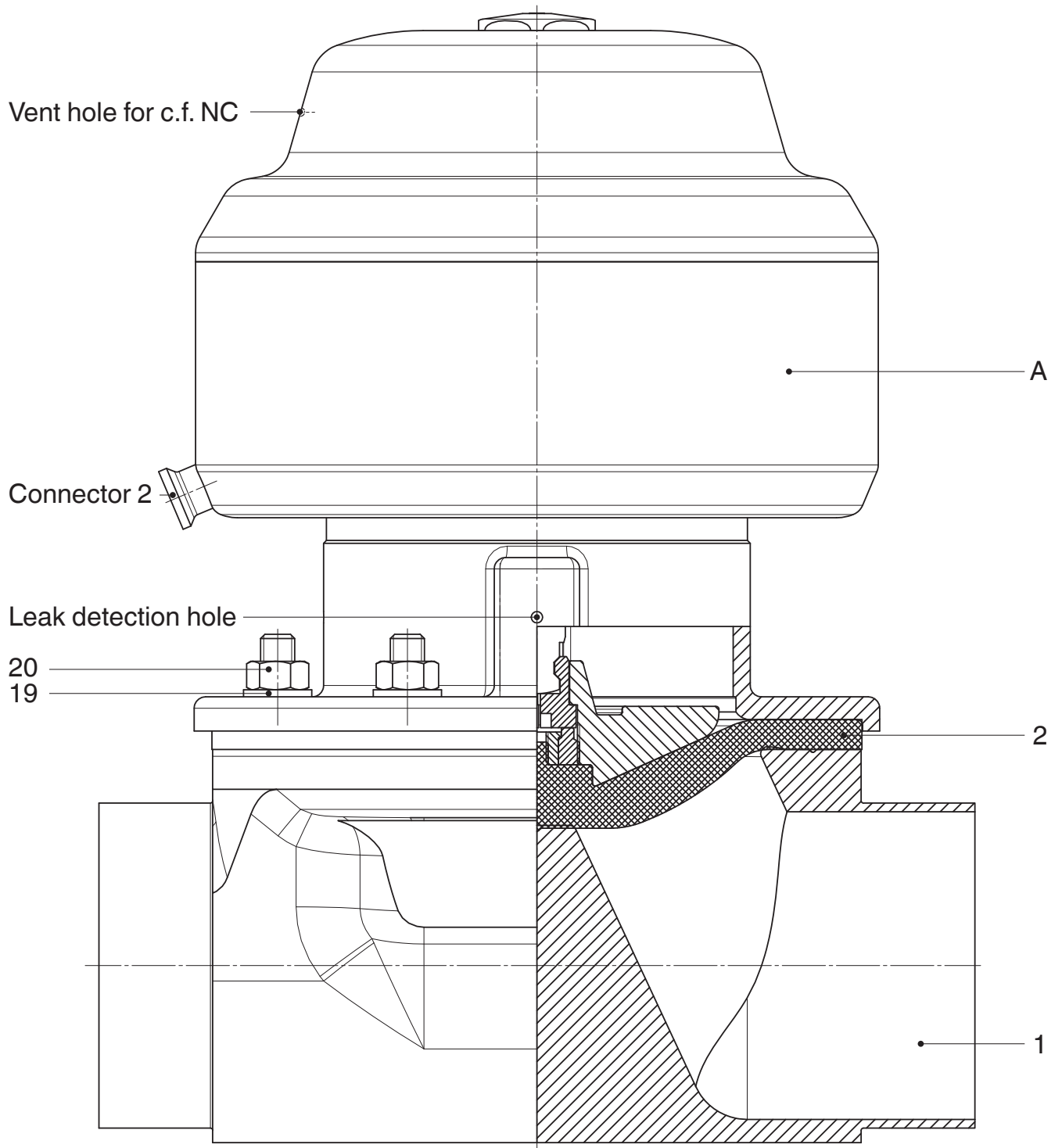Item	Name	Order description
Diaphragm size 100		
1	Valve body	K600...
2	Diaphragm	600...M
19	Washer	} 650...S30...
20	Nut	
A	Actuator	9650...

Declaration of Incorporation

according to the **EC Machinery Directive 2006/42/EC, Annex II, 1.B**
for partly completed machinery

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Description and identification of the partly completed machinery:

Make: GEMÜ Diaphragm valve, pneumatically operated
Serial number: from December 29, 2009
Project number: MV-Pneum-2009-12
Commercial name: Type 650

We hereby declare that the following essential requirements of the Machinery Directive 2006/42/EC have been fulfilled:

1.1.3., 1.1.5., 1.2.1., 1.3., 1.3.2., 1.3.3., 1.3.4., 1.3.7., 1.3.9., 1.5.3., 1.5.5., 1.5.6., 1.5.7., 1.5.8., 1.5.9., 1.6.5.

We also declare that the specific technical documentation has been compiled in accordance with part B of Annex VII.

We expressly declare that the partly completed machinery complies with the relevant provisions of the following EC directives:


2006/42/EC:2006-05-17: (Machinery Directive) Directive 2006/42/EC of the European Parliament and of the Council of 17 May 2006 on machinery, and amending Directive 95/16/EC (recast) (1)

The manufacturer or his authorised representative undertake to transmit, in response to a reasoned request by the national authorities, relevant information on the partly completed machinery. This transmission takes place:

electronically

This does not affect the intellectual property rights!

Important note! The partly completed machinery may be put into service only if it was determined, where appropriate, that the machinery into which the partly completed machinery is to be installed meets the provisions of this Directive.

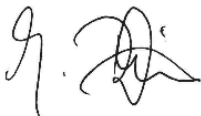Joachim Brien
Head of Technical Department

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declare that the equipment listed below complies with the safety requirements of the Pressure Equipment Directive 2014/68/EU.

Description of the equipment - product type

Diaphragm Valve
GEMÜ 650


Notified body: TÜV Rheinland
Berlin Brandenburg
Number: 0035
Certificate no.: 01 202 926/Q-02 0036
Applied standards: AD 2000

Conformity assessment procedure:
Module H

Note for equipment with a nominal size \leq DN 25:

The products are developed and produced according to GEMÜ process instructions and quality standards which comply with the requirements of ISO 9001 and of ISO 14001.

According to section 4, paragraph 3 of the Pressure Equipment Directive 2014/68/EU these products must not be identified by a CE-label.

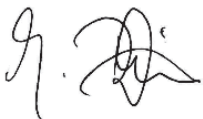Joachim Brien
Head of Technical Department

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Hereby we, **GEMÜ Gebr. Müller Apparatebau GmbH & Co. KG**